



ORagene[®]·DISCOVER


For collection of human DNA

REF OGR-500


For Research Use Only

Not for use in diagnostic procedures


1

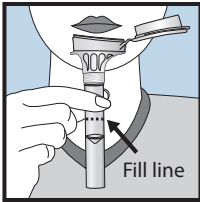


2

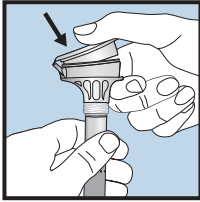


3

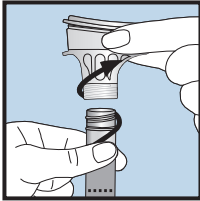

4

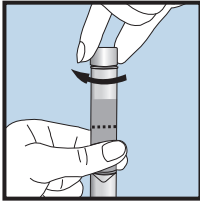

5

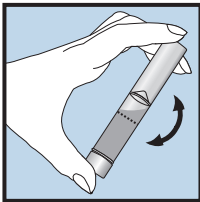

DNAgenotek




Made in Canada
DNA Genotek Inc.
Ottawa, ON, Canada K2K 1L1
Tel.: 613.723.5757
info@dnagenotek.com
www.dnagenotek.com

*Superior samples
Proven performance*

PD-PR-00157-ML Issue 4/2011-09

ENGLISH User instructions

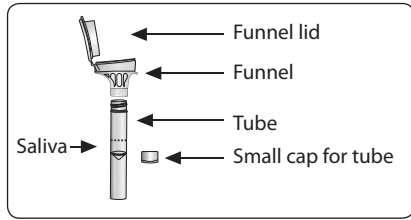

Read all instructions prior to collection.

This product is designed for the collection of human DNA from saliva samples.

Contents: Kit contains stabilizing liquid.

Warnings and precautions: Wash with water if stabilizing liquid comes in contact with eyes or skin. Do not ingest. See MSDS at www.dnagenotek.com.

Small cap, choking hazard.

Storage: 15°C / 30°C

Summary and explanation of the kit:

Oragene·DISCOVER is a self-collection kit that provides the materials and instructions for collecting and stabilizing saliva specimens.

Collection precautions:

Do NOT eat, drink, smoke or chew gum for 30 minutes before giving your saliva sample.

Do NOT remove the plastic film from the funnel lid.

Procedure:

Most people take between 2 and 5 minutes to deliver a saliva sample following steps 1 to 5.

- 1 Spit into funnel until the amount of liquid saliva (not bubbles) reaches the fill line shown in picture #1.
- 2 Hold the tube upright with one hand. Close the funnel lid with the other hand (as shown) by firmly pushing the lid until you hear a loud click. The liquid in the lid will be released into the tube to mix with the saliva. Make sure that the lid is closed tightly.
- 3 Hold the tube upright. Unscrew the funnel from the tube.
- 4 Use the small cap to close the tube tightly.
- 5 Shake the capped tube for 5 seconds. Discard or recycle the funnel.

Label legend:

- Consult package insert
- Collect saliva by (Use by)
- Catalog number
- Industrial Design Patent
- Caution, consult instructions for use
- Storage instructions
- Manufacturer